

Air-Dryable™ qPCR Mix

Product Handling Guide

Shipping:	On Dry/Blue Ice
Catalog number:	MDX082
Batch No.:	See vial
Concentration:	4x

Store at -20 °C



Storage and stability:

Air-Dryable™ qPCR Mix is shipped on dry/blue ice. On arrival store at -20 °C for optimum stability. Repeated freeze/thaw cycles should be avoided. Thawing during transportation does not affect the product performance. Solutions should be mixed/equilibrated after each thawing to avoid phasing.

Expiry:

When stored under the recommended conditions and handled correctly, full activity of the kit is retained until the expiry date on the outer box label.

Safety precautions:

Read and understand the SDS (Safety Data Sheets) before handling the reagents. Hardcopies of the SDS will be provided with the first shipment, thereafter they will be available upon request.

Quality control:

Meridian operates under ISO 13485 Quality Management System. Air-Dryable™ qPCR Mix and its components are extensively tested for activity, processivity, efficiency, heat activation sensitivity, absence of nuclease contamination and absence of nucleic acid contamination.

Notes:

For research or further manufactured use only.

Description

Air-Dryable™ qPCR Mix is a glycerol-free qPCR mix containing an antibody-mediated hot-start Taq polymerase, reaction buffer, dNTPs, MgCl₂ and air-dry compatible excipients. In order to produce room temperature air-dried qPCR reagents, assay specific primers and probes can be added to Air-Dryable™ qPCR Mix for subsequent air-drying.

Kit components

Table 1

Component
Air-Dryable™ qPCR Mix, 4x

Users Guidelines

Master mix preparation

Recommended reagent volumes of Air-dryable qPCR Mix and Primer-Probe Mix for air-drying are given in Table 2. Volumes are indicated per 20 µL final rehydrated reaction.

Table 2

Reagent	Volume
Air-Dryable™ qPCR Mix, 4x	5 µL
Primer-Probe Mix, 20x	1 µL

Dispense into reaction vessels, immediately transfer into a convection oven and run a suitable drying cycle.

For long-term storage at ambient temperatures, the air-dried product from the oven should be packaged with silica desiccant in a heat sealed pouch.

Assay setup

Rehydrate the air-dried qPCR master mix with 20 µL template-containing solution, vortex and run qPCR.

The qPCR conditions in Table 3 are suitable for amplicons of up to 200 bp. These cycling parameters are compatible with Air-Dryable™ qPCR Mix on a number of platforms, however they can be varied to suit different detection assay and machine-specific protocols.

Table 3

Step	Temperature	Time	Cycles
Polymerase activation	95 °C	2 min	1
Denaturation	95 °C	5 s	45
Annealing/Extension	60 °C	20 s	

Associated Products

Component	Cat. No.
Inhibitor-Tolerant qPCR Mix	MDX013
Air-Dryable™ Direct DNA qPCR Blood	MDX092
Air-Dryable™ Direct DNA qPCR Saliva	MDX130
Air-Dryable™ 1-Step RT-qPCR Mix	MDX095
Air-Dryable™ Direct DNA qPCR Plant	MDX116
Lyo-Ready qPCR Mix	MDX021
Lyo-Ready 1-Step RT-qPCR Mix	MDX024

Technical Support

For any technical enquiries, please contact our Technical Support team via email at: mbi.tech@meridianlifescience.com