

# Lyo-Ready™ Genotyping Direct qPCR Stool, 4x

## Product Handling Guide

Shipping:	Blue Ice
Catalog number:	MDX148
Batch No.:	See vial
Concentration:	4x

Store at -20 °C



### Storage and stability:

Lyo-Ready™ Genotyping Direct qPCR Stool is shipped on blue ice. On arrival store at -20 °C for optimum stability. Repeated freeze/thaw cycles should be avoided. Thawing during transportation does not affect the product performance. Solutions should be mixed/equilibrated after each thawing to avoid phasing.

### Expiry:

When stored under the recommended conditions and handled correctly, full activity of the kit is retained until the expiry date on the outer box label.

### Safety precautions:

Read and understand the SDS (Safety Data Sheets) before handling the reagents. Hardcopies of the SDS will be provided with the first shipment, thereafter they will be available upon request.

### Quality control:

Meridian operates under ISO 13485 Quality Management System. Lyo-Ready™ Genotyping Direct qPCR Stool and its components are extensively tested for activity, efficiency, sensitivity, absence of nuclease contamination and absence of nucleic acid contamination.

### Notes:

For research or further manufactured use only.

## Description

Lyo-Ready™ Genotyping Direct qPCR Stool is a glycerol-free, one tube formulation compatible with all dual-label probe and intercalating dye chemistries for detection of genetic variants, such as single nucleotide variants and copy number variants. The formulation combines the latest advances in buffer chemistry and PCR enhancers, together with an optimized concentration of antibody-mediated hot-start polymerase, dNTPs and MgCl<sub>2</sub>. Lyo-Ready™ Genotyping Direct qPCR Stool has been designed for fast, precise, and highly reproducible allelic discrimination and cluster separation with SNP detection assays, utilising allele specific probe and/or primer methods, even in the presence of PCR inhibitors like crude Stool sample. In order to produce room temperature lyophilized qPCR reagents, assay specific primers and probes can be added to Lyo-Ready™ Genotyping Direct qPCR Stool for subsequent lyophilization.

## Kit components

Table 1

Component
Lyo-Ready™ Genotyping Direct qPCR Stool, 4x

## Users Guidelines

### Inhibitor Tolerance

The amount of inhibition tolerated by Lyo-Ready™ Genotyping Direct qPCR Stool is variable depending on several factors, including assay design and sample quality. For this reason, an initial sample titration is recommended.

Tolerance can be further improved by heating stool in water or buffer at 95°C for 15 minutes before adding to reactions.

### Master mix preparation

Recommended reagent volumes of Lyo-Ready™ Genotyping Direct qPCR Stool and Primer–Probe Mix for lyophilization are given in Table 2. Volumes are indicated per 20 µL final rehydrated reaction.

Table 2

Reagent	Volume
Lyo-Ready™ Genotyping Direct qPCR Stool, 4x	5 µL
Primer–Probe Mix, 20x	1 µL*
Water	x µL
<b>Total Volume</b>	<b>Up to 20 µL</b>

\*Primer and probe concentration needs to be optimised

Dispense into reaction vessels, immediately transfer into a freeze-drier and run a suitable drying cycle.

### Lyophilization

For lyophilization protocols, please consult our “Lyophilization & Post-Lyophilization User Guideline”.

For long-term storage at ambient temperatures, the lyophilized product from the freeze-dryer should be packaged with a silica sachet in a heat sealed pouch at low relative humidity conditions.

## Assay setup

Rehydrate the lyophilized qPCR master mix in the reaction vials with 20 µL template-containing solution and run qPCR.

The qPCR conditions in Table 3 are suitable for amplicons of up to 200 bp. These cycling parameters have been optimized for Lyo-Ready™ Genotyping Direct qPCR Stool on a number of platforms, however they can be varied to suit different machine-specific protocols.

Table 3

Step	Temperature	Time	Cycles
Polymerase activation	95 °C	5 min	1
Denaturation	95 °C	15 s	45
Annealing/Extension**	60-68 °C	30 s	

\*\*Annealing/Extension temperature needs to be optimised based on probes used.

## Associated Products

Product	Cat. No.
Lyo-Ready™ Genotyping Direct qPCR Blood, 4x	MDX128
Lyo-Ready™ Genotyping Direct qPCR Urine, 4x	MDX158
Lyo-Ready™ Genotyping Direct qPCR FFPE, 4x	MDX168
Air-Dryable™ Direct DNA qPCR Stool	MDX140
Air-Dryable™ Direct RNA/DNA qPCR Stool	MDX141
Lyo-Ready™ Direct DNA qPCR Stool	MDX142
Lyo-Ready™ Direct RNA/DNA qPCR Stool	MDX143

## Technical Support

For any technical enquiries, please contact our Technical Support team via email at: [mbi.tech@meridianlifescience.com](mailto:mbi.tech@meridianlifescience.com)